

CHANGING PROCEDURES FOR STERILE CLEANROOMS

Make sure that you have the correct garment specification and fabric for your manufacturing environment
All Cleanroom clothing should be processed to the correct standard in a dedicated Cleanroom laundry facility

<p>1</p>  <p>NO Cosmetics NO Jewellery NO Watches NO Personal Items</p>	<p>2</p>  <p>All operators should wear a polyester undergarment</p>	<p>3</p>  <p>Put on a mob cap and shoe covers</p>	<p>4</p>  <p>Wash hands</p>
<p>5</p>  <p>Put on first pair of sterile gloves</p>	<p>6</p>  <p>Select cleanroom garments</p>	<p>7</p>  <p>Put on face mask ensuring a good fit</p>	<p>8</p>  <p>Put on hood</p>
<p>9</p>  <p>Hold coverall at waist level making sure it does not touch the floor</p>	<p>10</p>  <p>Put coverall on and tuck the hood in</p>	<p>11</p>  <p>Sit on the step over bench, put on the first overboot and swing leg to the clean side of the bench. Repeat with second overboot</p>	<p>12</p>  <p>Put on goggles</p>
<p>13</p>  <p>Put on second pair of sterile gloves</p>	<p>14</p>  <p>Check in the mirror and enter the cleanroom</p>	<p>15</p>  <p>To disrobe please follow the steps in reverse order</p>	<p>16</p>  <p>All used garments should be isolated and sent to a Cleanroom laundry facility for processing</p>

info@cmsupply.co.uk

www.cmsupply.co.uk